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## **Safety instructions (E)**

### **1) Work Protection and Safety**

When handling and working with Dewar containers, the use of safety goggles and protective gloves is necessary. Regional specific safety regulations referring to glass apparatuses, liquefied gasses, laboratory devices, vacuum insulated glass containers, etc., must be strictly followed.

### **2) Regulation and Standards**

KGW-ISOTHERM Dewar containers meet the standards of DIN 12492.

They are manufactured in consideration of the European Pressure Specifications 97/23 EG from Borosilicate glass 3.3 under DIN/ISO 3585 DURAN (AD-data sheet N4).

Standard Dewar containers are allowed to be used up to just under 0.1 bar pressure (Inquiries about higher pressure requirements are welcome).

Dewar vessels with tight-closable lid have to be provided with a gas-release hole of a 2mm diameter, if liquid gas or dry ice is used. Dewar containers made of glass are to be used with their protective jackets only (example: metal safety cover). Unjacketed Dewar containers (Type A), are only allowed in the internal construction of an apparatus or as a replacement piece. Safety glasses and protective gloves are a necessity when installing Type A Dewar container in a metal jacket or in an apparatus.

### **3) Inspection before Use**

Before using a Dewar container, a complete inspection must be made for scratches, cracks, chips or other flaws. Containers with surface flaws that will be subjected to mechanical or thermal stresses cannot be used due to the threat of implosion.

### **4) Prepare for Use**

To preserve the function, quality and safety of a Dewar container, special attention should be paid that the container is used only in the temperature range of -200°C to +150°C. Dewar containers that are to be used beyond these temperature limits must undergo additional manufacturing processes.

Standard Dewar containers of glass cannot be used to contain liquid or gaseous Helium. Higher technical specifications are required for the containment of Helium and are available upon request.

### **5) Maintenance and Cleaning**

Dewar containers are maintenance-free. By the eventual cleaning of a Dewar glass container, special care must be taken not to damage the surface. The glass container cannot be cleaned with metal or ceramic objects.

We are always available for our customers to answer any questions.

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